

THE INSTITUTE FOR GENOMIC RESEARCH
Standard Operating Procedure

TITLE: **AMINOALLYL LABELING OF RNA FOR MICROARRAYS**

PAGE: 1 of 7

SOP #: M004

REVISION LEVEL: 1

EFFECTIVE DATE: 6/21/01

AUTHOR:

Jeremy Hasseman

PRIMARY REVIEWERS:

Emily Chen, Erik Snesrud, Ivana Yang

1. PURPOSE

This protocol describes the labeling of eukaryotic RNA with aminoallyl labeled nucleotides via first strand cDNA synthesis followed by a coupling of the aminoallyl groups to either Cyanine 3 or 5 (Cy 3/Cy5) fluorescent molecules.

2. SCOPE

This procedural format is utilized by Human Colon Cancer and Mouse microarray projects under the supervision of John Quackenbush within the Eukaryotic Genomics Dept.

3. MATERIALS

- 3.1 5-(3-aminoallyl)-2'-deoxyuridine-5'-triphosphate (AA-dUTP) (Sigma; Cat # A0410)
- 3.2 100 mM dNTP Set PCR grade (Life Technologies; Cat # 10297-018)
- 3.3 Random Hexamer primers (3mg/mL) (Life Technologies; Cat # 48190-011)
- 3.4 SuperScript II RT (200U/μL) (Life Technologies; Cat # 18064-014)
- 3.5 Cy-3 ester (AmershamPharmacia; Cat # PA23001)
- 3.6 Cy-5 ester (AmershamPharmacia; Cat # PA25001)
- 3.7 QIAquick PCR Purification Kit (Qiagen; Cat # 28106)
- 3.8 RNeasy[®] Mini Kit (Qiagen; Cat # 74106)

4. REAGENT PREPARATION

4.1 Phosphate Buffers

4.1.1 Prepare 2 solutions: 1M K₂HPO₄ and 1M KH₂PO₄

4.1.2 To make a 1M Phosphate buffer (KPO₄, pH 8.5-8.7) combine:

1M K₂HPO₄.....9.5 mL

1M KH₂PO₄.....0.5 mL

THE INSTITUTE FOR GENOMIC RESEARCH

TITLE: AMINOALLYL LABELING OF RNA FOR MICROARRAYS

SOP #: M004

REVISION LEVEL: 0

PAGE: 2 of 7

- 4.1.3 For 100 mL Phosphate wash buffer (5 mM KPO₄, pH 8.0, 80% EtOH) mix:

1 M KPO₄ pH 8.5..... 0.5 mL
 MilliQ water.....15.25 mL
 95% ethanol..... 84.25 mL

Note: Wash buffer will be slightly cloudy.

- 4.1.4 Phosphate elution buffer is made by diluting 1 M KPO₄, pH 8.5 to 4 mM with MilliQ water.

4.2 Aminoallyl dUTP

- 4.2.1 For a final concentration of 100mM add 19.1 µL of 0.1 M KPO₄ buffer (pH 7.5) to a stock vial containing 1 mg of aa-dUTP. Gently vortex to mix and transfer the aa-dUTP solution into a new microfuge tube. Store at –20°C.

- 4.2.2 Measure the concentration of the aa-dUTP solution by diluting an aliquot 1:5000 in 0.1 M KPO₄ (pH 7.5) and measuring the OD₂₈₉. (Stock concentration in mM = OD₂₈₉ x 704)

4.3 Labeling Mix (50X) with 2:3 aa-dUTP: dTTP ratio

- 4.3.1 Mix the following reagents:

	Final concentration
dATP (100 mM).....	5µL.....(25 mM)
dCTP (100 mM).....	5µL.....(25 mM)
dGTP (100 mM).....	5µL.....(25 mM)
dTTP (100 mM).....	3µL.....(15 mM)
aa-dUTP (100 mM).....	2µL.....(10 mM)
Total:	20µL

- 4.3.1 Store unused solution at –20°C.

4.4 Sodium Carbonate Buffer (Na₂CO₃): 1M, pH 9.0

- 4.4.1 Dissolve 10.8 g Na₂CO₃ in 80 mL of MilliQ water and adjust pH to 9.0 with 12 N HCl; bring volume up to 100 mL with MilliQ water.

TITLE: AMINOALLYL LABELING OF RNA FOR MICROARRAYS

SOP #: M004

REVISION LEVEL: 0

PAGE: 3 of 7

4.4.2 To make a 0.1 M solution for the dye coupling reaction dilute 1:10 with water.

Note: Carbonate buffer changes composition over time; make it fresh every couple of weeks to a month.

4.5 Cy-dye esters

4.5.1 Cy3-ester and Cy5-ester are provided as a dried product in 5 tubes. Resuspend a tube of dye ester in 73 μ L of DMSO before use.

4.5.2 Wrap all reaction tubes with foil and keep covered as much as possible in order to prevent photobleaching of the dyes.

Note: Dye esters must either be used immediately or aliquoted and stored at -80°C . Any introduced water to the dye esters will result in a lower coupling efficiency due to the hydrolysis of the dye esters. Since DMSO is hygroscopic (absorbs water from the atmosphere) store it well sealed in desiccant.

5. PROCEDURE

5.1 Aminoallyl Labeling

5.1.1 To 10 μ g of total RNA (or 2 μ g poly(A⁺) RNA) which has been DNase I-treated and Qiagen RNeasy purified, add 2 μ L Random Hexamer primers (3mg/mL) and bring the final volume up to 18.5 μ L with RNase-free water.

5.1.2 Mix well and incubate at 70°C for 10 minutes.

5.1.3 Snap-freeze in dry ice/ethanol bath for 30 seconds, centrifuge briefly at $>10,000$ rpm and continue at room temperature.

5.1.4 Add:

5X First Strand buffer..... 6 μ L
0.1 M DTT..... 3 μ L
50X aminoallyl-dNTP mix..... 0.6 μ L
SuperScript II RT (200U/ μ L)..... 2 μ L

TITLE: AMINOALLYL LABELING OF RNA FOR MICROARRAYS

SOP #: M004

REVISION LEVEL: 0

PAGE: 4 of 7

5.1.5 Mix and incubate at 42°C for 3 hours to overnight.

5.1.6 To hydrolyze RNA, add:

1 M NaOH 10 µL
0.5 M EDTA 10 µL

mix and incubate at 65°C for 15 minutes.

5.1.7 Add 10 µL of 1 M HCl to neutralize pH. (Alternatively, one can add 25 µL 1 M HEPES pH 7.0 or 25 µL 1 M Tris pH 7.4)

5.2 Reaction Purification I: Removal of unincorporated aa-dUTP and free amines (Qiagen PCR Purification Kit)

Note: This purification protocol is modified from the Qiagen QIAquick PCR purification kit protocol. The phosphate wash and elution buffers (prepared in 4.1.3 & 4.1.4) are substituted for the Qiagen supplied buffers because the Qiagen buffers contain free amines which compete with the Cy dye coupling reaction.

5.2.1 Mix cDNA reaction with 300 µL (5X reaction volume) buffer PB (Qiagen supplied) and transfer to QIAquick column.

5.2.2 Place the column in a 2 ml collection tube (Qiagen supplied) and centrifuge at ~13,000 rpm for 1 minute. Empty collection tube.

5.2.3 To wash, add 750 µL phosphate wash buffer to the column and centrifuge at ~13,000 rpm for 1 minute.

5.2.4 Empty the collection tube and repeat the wash and centrifugation step (5.2.3).

5.2.5 Empty the collection tube and centrifuge column an additional 1 minute at maximum speed.

5.2.6 Transfer column to a new 1.5 mL microfuge tube and carefully add 30 µL phosphate elution buffer (*see 4.1.4*) to the center of the column membrane.

TITLE: AMINOALLYL LABELING OF RNA FOR MICROARRAYS

SOP #: M004

REVISION LEVEL: 0

PAGE: 5 of 7

5.2.7 Incubate for 1 minute at room temperature.

5.2.8 Elute by centrifugation at ~13,000 rpm for 1 minute.

5.2.9 Elute a second time into the same tube by repeating steps 5.2.6-5.2.8. The final elution volume should be ~60 µL.

5.2.10 Dry sample in a speed vac.

5.3 Coupling aa-cDNA to Cy Dye Ester.

5.3.1 Resuspend aminoallyl-labeled cDNA in 4.5 µL 0.1 M sodium carbonate buffer (Na₂CO₃), pH 9.0.

Note: Carbonate buffer changes composition over time so make sure you make it fresh every couple of weeks to a month.

5.3.2 Add 4.5 µL of the appropriate NHS-ester Cy dye (prepared in DMSO: *see 4.5*)

Note: To prevent photobleaching of the Cy dyes wrap all reaction tubes in foil and keep them sequestered from light as much as possible.

5.3.3 Incubate the reaction for 1 hour in the dark at room temperature.

5.4 Reaction Purification II: Removal of uncoupled dye (Qiagen PCR Purification Kit)

5.4.1 To the reaction add 35 µL 100 mM NaOAc pH 5.2.

5.4.2 Add 250 µL (5X reaction volume) Buffer PB (Qiagen supplied).

5.4.3 Place a QIAquick spin column in a 2 mL collection tube (Qiagen supplied), apply the sample to the column, and centrifuge at ~13,000 for 1 minute. Empty collection tube.

5.4.4 To wash, add 0.75 mL Buffer PE (Qiagen supplied) to the column and centrifuge at ~13,000 for 1 minute.

Note: Make sure Buffer PE has added ethanol before using (see label for correct volume).

TITLE: AMINOALLYL LABELING OF RNA FOR MICROARRAYS

SOP #: M004

REVISION LEVEL: 0

PAGE: 6 of 7

- 5.4.5 Empty collection tube and centrifuge column for an additional 1 minute at maximum speed.
- 5.4.6 Place column in a clean 1.5 mL microfuge tube and carefully add 30 µL Buffer EB (Qiagen supplied) to the center of the column membrane.
- 5.4.7 Incubate for 1 minute at room temperature.
- 5.4.8 Elute by centrifugation at ~13,000 rpm for 1 minute.
- 5.4.9 Elute a second time into the same tube by repeating steps 5.4.6-5.4.8. The final elution volume should be ~60 µL.

Note: This protocol is modified from the Qiagen QIAquick Spin Handbook (04/2000, pg. 18).

5.5 Analysis of Labeling Reaction

- 5.5.1 Use a 50 µL Beckman quartz MicroCuvette to analyze the entire undiluted sample in a spectrophotometer.
- 5.5.2 Wash the cuvette with water and blow dry with compressed air duster.
- 5.5.3 Pipette sample into cuvette and place cuvette in spectrophotometer.
- 5.5.4 For each sample measure absorbance at 260 nm and either 550 nm for Cy3 or 650 nm for Cy5, as appropriate.
- 5.5.5 Pipette sample from cuvette back into the original sample tube.
- 5.5.6 For each sample calculate the total picomoles of cDNA synthesized using:

$$\text{pmol nucleotides} = \frac{[\text{OD}_{260} * \text{volume } (\mu\text{L}) * 37 \text{ ng}/\mu\text{L} * 1000 \text{ pg}/\text{ng}]}{324.5 \text{ pg}/\text{pmol}}$$

Note: 1 OD₂₆₀ = 37 ng/µL for cDNA; 324.5 pg/pmol average molecular weight of a dNTP)

TITLE: AMINOALLYL LABELING OF RNA FOR MICROARRAYS

SOP #: M004

REVISION LEVEL: 0

PAGE: 7 of 7

- 5.5.7 For each sample calculate the total picomoles of dye incorporation (Cy3 or Cy5 accordingly) using:

$$\text{pmol Cy3} = \frac{\text{OD}_{550} * \text{volume } (\mu\text{L})}{0.15}$$

$$\text{pmol Cy5} = \frac{\text{OD}_{650} * \text{volume } (\mu\text{L})}{0.25}$$

$$\text{nucleotides/dye ratio} = \frac{\text{pmol cDNA}}{\text{pmol Cy dye}}$$

Note: >200 pmol of dye incorporation per sample and a ratio of less than 50 nucleotides/dye molecules is optimal for hybridizations (see Microarray Cookbook II)

- 5.5.8 After analysis mix together the two differentially labeled probes (Cy3 vs. Cy5) which will be hybridized to the same microarray slide for study of relative gene expression.
- 5.5.9 Dry the Cy3/Cy5 probe mixture to completion in a speed vac and continue with SOP: M005 for the hybridization of the probe to a microarray.